

## BOTTLE-TOP DISPENSER PROBLEM DETERMINATION

<b>Problem</b>	<b>Probable cause</b>	<b>Chek point</b>
No suction of liquid	Is the inlet tube connection loose?	Tighten.
	Has the ball of the inlet valve or outlet valve become stuck?	Remove the inlet valve or outlet valve and cleaning.(See the page 9)
	Is there dirt on the inlet valve or outlet valve?	
Air bubbles	Is the inlet tube connector loose?	Tighten.
	Is the outlet valve loose?	Tighten with the provided wrench.
Volume indicator shift	Is the set screw loose?	Tighten.
Liquid leakage from the outlet valve.	Is the outlet valve loose?	Tighten with the provided wrench.
	Is the dispensing nozzle holder loose?	Tighten.
No discharge of liquid	Has crystallized liquid adhered to the inlet valve or outlet valve?	Cleaning. (See the page 9)

If the problem persists even after the above points have been checked, stop usage of the unit immediately and ask the distributor to repair it.

At this time please confirm that there has been no contamination by chemicals that are harmful to the body.